

20000016465

LR 1000.62 Adapter_052017

LR 1000.62 Adapter



Beachten Sie die Betriebsanleitung des LR 1000 basic/control!
Observe the Operating instructions of the LR 1000 basic/control!

Ursprungssprache

DE

Sicherheitshinweise

- Der Adapter **LR 1000.62** ist nur für den Deckel der **IKA LR 1000** Reaktorgefäße geeignet.
- Achten Sie auf den richtigen Sitz des Deckels in der Halterung. Prüfen Sie dies in regelmäßigen Abständen.
- Beachten Sie die Betriebsanleitung der Temperaturmessfühler.

- Verwenden Sie nur original **IKA** Ersatz- und Zubehörteile!
- Achten Sie auf einen korrekten Sitz der O-Ringe. Die O-Ringe dürfen nicht gequetscht oder beschädigt sein.

Zugelassene IKA Temperaturmessfühler:

- IKA Thermometer (Ø6 mm)

Source language: German

EN

Safety instructions


- The adapter **LR 1000.70** is only suitable for the lid of the reactor vessels of **IKA LR 1000**.
- Ensure correct fit of the lid in the holder. Check the tight fit at regular intervals.
- Refer to the operating instructions for the temperature sensors.

- Use only original **IKA** spare parts and accessories!
- Observe that the O-rings fit properly and aren't damaged or squeezed.

Permitted IKA temperature sensors:

- IKA Thermometer (Ø6 mm)

Montage / Assembling

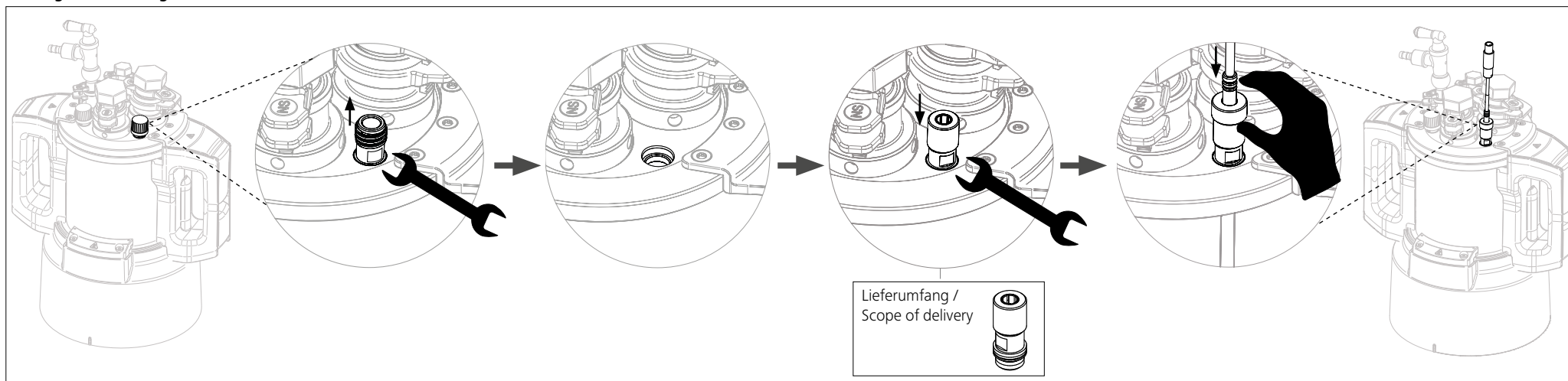

IKA®-Werke GmbH & Co.KG

Janke & Kunkel-Str. 10


D-79219 Staufen

Tel. +49 7633 831-0

Fax +49 7633 831-98

sales@ika.de

www.ika.com

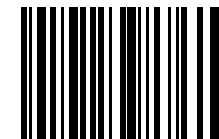

20014403