designed for scientists

ROTASTAND

5

/// ENGLISH

13

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21

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9

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17

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25

/// INDONESIA

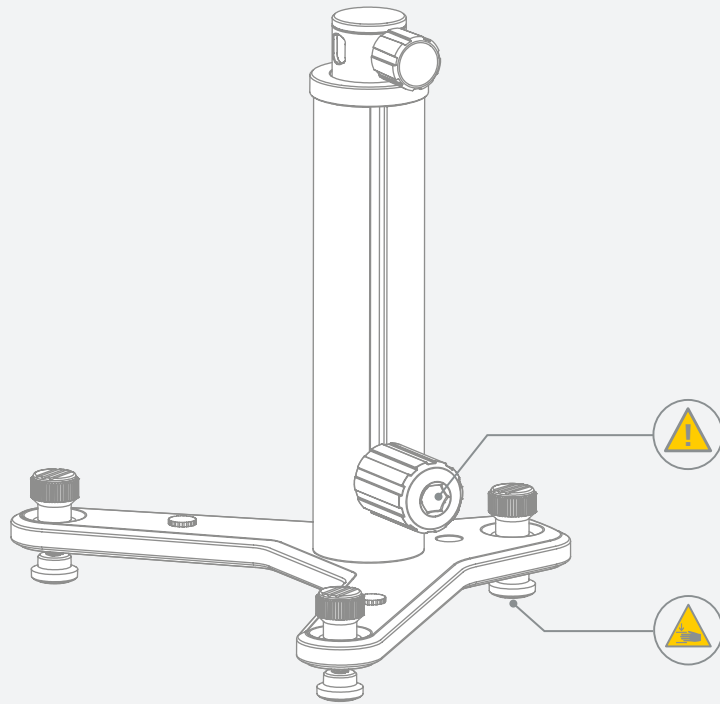
Fig. 1

	Explication of warning symbols	6
	Safety instructions.....	6
	Intended use.....	7
	Unpacking	7
	Assembly	7
	Maintenance and cleaning.....	8
	Technical Data	8
	Warranty	8

Source language: German

Explication of warning symbols

/// Warning symbols

Notice!

Indicates practices which, if not avoided, can result in equipment damage.

Caution!

Indicates crushing risk to fingers / hands.

/// General Symbols

A — Position number

Indicates device components relevant to actions.

Safety instructions

/// General information

- › **Read the operating instructions completely before starting up and follow the safety instructions. If this accessory is used with another device, observe also its operating instructions.**
- › Keep the operating instructions in a place where it can be accessed by everyone.
- › Follow the safety instructions, guidelines, occupational health, safety and accident prevention regulations.

Notice!

- › Pay attention to the marked sites in **Fig. 1**.

/// Device setup

- › Set up the device in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.
- › All screw connections must be properly tightened.
- › Check the mounting position of the viscometer to avoid the danger of the unit tipping over.

Caution!

- › Ensure that your hands / fingers do not get crushed when setting down the device.

/// Working with the device

Notice!

- › Risk of tipping exists when tilted the stand at an angle of being over 10° with full load.
- › Observe the maximum permissible load-bearing capacity (see "Technical data").
- › When moving the stand, ensure the device is at the lowest end stop position.
- › When used as intended and in accordance with the operating instructions, there is no danger to people, the environment or the device. The manufacturer cannot be held responsible for any damages as a result of incorrect use of the device or non-compliance with the safety instructions.

/// Accessories

- › Protect the device and accessories from bumping and impacting.
- › Check the device and accessories beforehand for damage each time when you use them. Do not use damaged components.

Intended use

/// Use

- › The IKA ROTASTAND has been designed for holding, lifting and lowering IKA ROTAVISC.

/// Area of use

- › Indoor environments similar to that a laboratory of research, teaching, trade or industry area.
- › The safety of the user cannot be guaranteed:
 - If the device is operated with accessories that are not supplied or recommended by the manufacturer.
 - If the device is operated improperly or contrary to the manufacturer's specifications.
 - If the device is modified by the third parties.


Unpacking

/// Unpacking



- › Unpack the device carefully. Any damage should immediately be reported to the carrier (mail, rail or freight forwarding company).

/// Scope of delivery

1	ROTASTAND pillar
2	ROTASTAND base
3	Screw washer
4	Hexagon socket screw
5	Hexagon socket screw (3x)
6	Screwdriver (AF = 2.5 mm) DIN ISO 2936
7	Screwdriver (AF = 5 mm) DIN ISO 2936

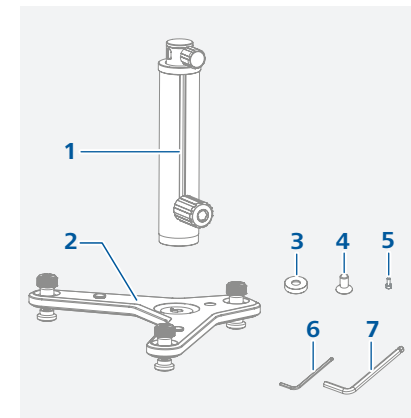Assembly

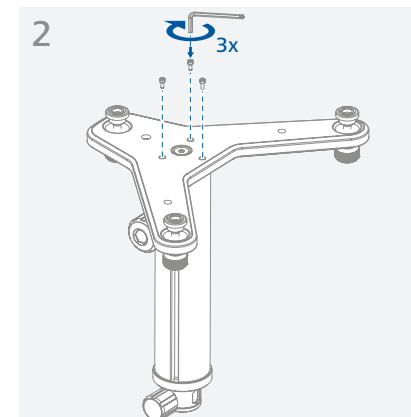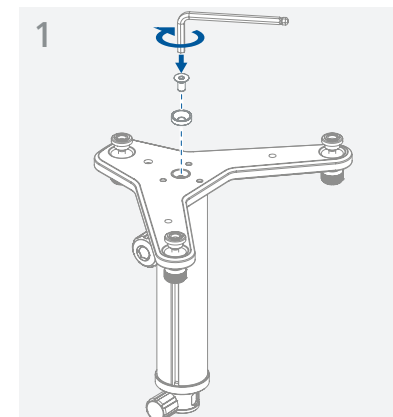Maintenance and cleaning

- › The device is maintenance-free. It is only subject to the natural wear and tear of components and their statistical failure rate.

/// Cleaning

- › Use only cleaning agents which have been approved by IKA to clean the devices:
These are: water (containing surfactant) and isopropyl alcohol.
- › Wear protective gloves during cleaning the devices.
- › Do not allow moisture to get into the device when cleaning.
- › Before using another than the recommended method for cleaning or decontamination, the user must ascertain with IKA that this method does not destroy the device.

/// Ordering spare parts

- › When ordering spare parts, please give:
 - Device type,
 - position number and description of spare part, see www.ika.com.

/// Repairs

- › Please only send devices in for repair that have been cleaned and are free of materials which might present health hazards.
- › For repair, please request the “**Safety Declaration (Decontamination Certificate)**” from IKA or use the download printout of it from IKA website at www.ika.com.
- › If your appliance requires repair, return it in its original packaging. Storage packaging is not sufficient when sending the device - also use appropriate transport packaging.

Technical data

Stroke max.	mm	200
Diameter (Ø)	mm	16
Dynamic load	kg	5
Dimensions (W x D x H)	mm	351 x 246 x 318
Weight	kg	4.8

› Subject to technical changes!

Warranty

- › In accordance with IKA Terms and Conditions of Sale, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.
- › The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

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IKA India Private Limited
IKA England LTD.
IKA Vietnam Company Limited
IKA Works (Thailand) Co. Ltd.